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WindowsDirect 電腦直傳

### Customers in the Food Product Industry

- Research and development of food products, drinks, and food additives.
- Inspections of food products manufacturing processes and products.
- Measurements of raw materials (such as rice, malt, tea leaves, and corn starch)

### Customers in the Pharmaceuticals and Cosmetics Industries

- Pharmaceuticals research and development (pills, granules, capsules, and ointments)
- Cosmetics research and development (such as hair products, face lotions, and soap)
- Inspections of pharmaceutical and cosmetic manufacturing processes and products

### Customers in Water Supply and Environmental Fields

- Moisture ratio measurements and inspections of sludge
- Moisture ratio measurements and inspections of soil
- Moisture ratio measurements and inspections of waste matter

### Customers in the Chemical Industry

- Research and development of samples, and inspection of products (solids, powders, pellets, films, and liquid samples)
- Moisture ratio measurements for plastics, rubbers, paints, pesticides, and functional materials

### Customers in Metals and Raw Materials Industries

- Moisture ratio measurements of raw materials (such as iron ore)
- Moisture ratio measurements of metals (such as metal powders)
- Moisture ratio measurements of lumber, paper, fiber, and lime

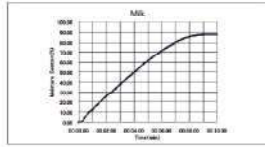
### Customers in the Automotive and Electrical Machinery Industries

- Moisture ratio measurements of plastic pellets and plastic parts
- Moisture ratio measurements of toner
- Moisture ratio measurements of foundry sand

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The drying curve for milk in timed ending mode is shown below.


Photos of the milk before and after drying are shown below.

(Before measurement) 1 g of milk was dripped on to a fibreglass sheet for liquid measurements.  
(After measurement) The moisture has evaporated from the milk, and the remaining fat has yellowed slightly.

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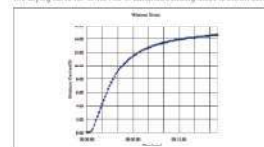The drying curve for white rice in automatic ending mode is shown below.


Photos of the white rice before and after drying are shown below.


(Before measurement) The white rice was spread evenly over the pan.  
(After measurement) The entire sample turned yellow.

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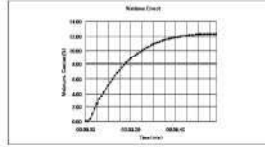The drying curve for corn starch in automatic ending mode is shown below.

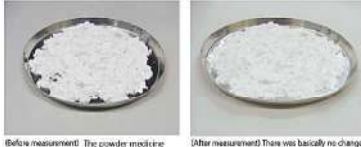
Powder medicine: before and after drying are shown below.

(Before measurement) The powder medicine was spread evenly over the pan.  
(After measurement) There was basically no change in appearance.

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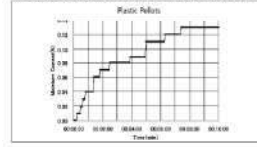The drying curve for plastic pellets in timed ending mode is shown below.

Photos of the plastic pellets before and after drying are shown below.

(Before measurement) The sample was spread evenly over the pan.  
(After measurement) There was basically no change in appearance.

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Sample	Sample Quantity	Measurement Mode		Set Temperature (°C)	Measurement Time (min)	Moisture Ratio (%)
		Ending Conditions	Finishing Conditions (% or min)			
Dog food	1g	AUTO	0.05%	160	5:48	6.45
Table salt	5g	TIME	10 minutes	200	10:00	0.08
Instant coffee	1g	TIME	10 minutes	120	10:00	7.43
Coffee beans (raw)	5g	AUTO	0.05%	140	17:30	9.32
Coffee beans (roasted)	3g	AUTO	0.05%	140	7:06	2.68
Green tea	5g	AUTO	0.05%	120	9:05	3.76
Corn starch	5g	AUTO	0.02%	180	9:25	12.17
Sugar (granulated sugar)	5g	AUTO	0.05%	160	1:02	0.13
White rice	6g	AUTO	0.05%	200	13:55	14.48
Mayonnaise	1g	TIME	10 minutes	160	10:00	20.51
Orange juice	1g	AUTO	0.05%	140	10:09	88.89
Milk	1g	AUTO	0.05%	140	7:30	87.36
Chocolate	3g	AUTO	0.01%	140	6:18	2.36
Rolled oats	6g	AUTO	0.05%	200	10:05	12.65
Tomato ketchup	2.5g	AUTO	0.1%	140	19:47	69.40
Frozen sweets	2.5g	TIME	12 minutes	140	12:00	84.53
Dried mangoes	5g	AUTO	0.05%	120	28:27	6.62
Palm oil	2.5g	TIME	5 minutes	120	5:00	0.41
Hand soap	1g	AUTO	0.05%	200	21:36	88.89
Lipstick	1g	TIME	3 minutes	100	3:00	0.73
Plastic (PMMA pellet)	10g	TIME	25 minutes	100	25:00	0.13
Photocopier paper	1g	AUTO	0.05%	200	1:50	7.84
Sodium tartrate dihydrate	5g	TIME	15 minutes	160	15:00	15.80
Detergent (powdered)	5g	AUTO	0.05%	160	13:08	9.79
Solid soap	3g	TIME	16 minutes	200	16:00	9.09
Water-based paint	1g	AUTO	0.05%	200	9:27	52.39
Sludge cake	2g	AUTO	0.05%	200	21:31	81.55
Potting soil	5g	AUTO	0.05%	120	15:30	33.40
Sawdust	4g	AUTO	0.05%	160	8:27	34.38